

IMPORTANT PLEASE READ BEFORE USING THIS COLUMN

Dear EZ Flash II User,

In order to improve serviceability and reduce the vast range of column types ACS (Advanced Chromatography Systems) and CSI (Cambridge Scientific Instruments) have changed the column winding to a new common column for all instruments.

The new column has longer arms and is to be mounted on the GC columns holder. The longer arms allow each end of the column to be assembled and installed separately.

This allows the user to check column flow on installation and gives the ability to change the fused silica tails more easily when they become contaminated.

See over for installation instructions. If any more help is required with the new column installation please contact us.

Advanced Chromatography Systems
3547 Meeks Farm Road
Suite D-1
Charleston
SC 29455
USA

Tel: 843 330 3538

Fax: 843 746 4545

Email: jam@advancedlcgcs.com

Website: <http://www.advancedlcgcs.com>



Cambridge Scientific Instruments
12-15 Sedgeway Business Park
Witchford
Cambridgeshire
CB6 2HY
UK

Tel: +44 (0)1353 669916

Fax: +44 (0)1353 669917

Email: camsci@btconnect.com


Website: <http://www.camsci.co.uk>

EZ Flash II Column Installation Guide

1



Attach the column to the column hanger as shown by fastening it with the wire supplied.

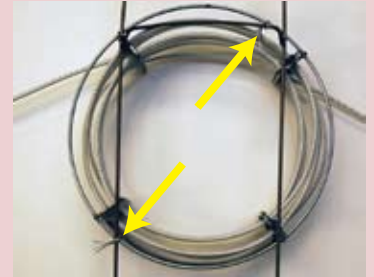

2



The column can now be hanged in the GC oven.


3

The end column end fittings can now be added.



4

Insert the end of the column extending from the right hand side round and underneath into the injector as shown.

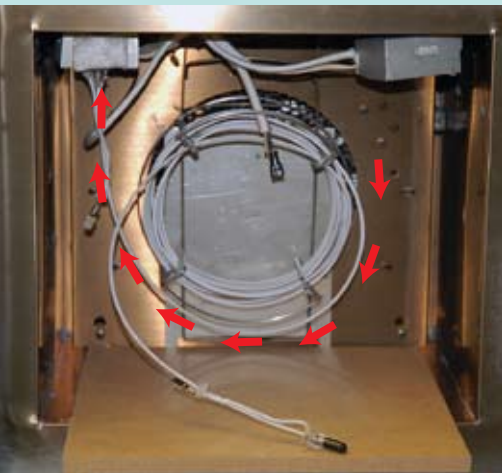

5

Finally prepare the other end fitting and insert the end of the column extending from the left hand side into the detector as shown.

